

ARCIS DNA Sample Prep (Bulk Kit)

UFL002 ARCIS DNA Sample Prep 48 rxn

Instructions for use

1. General Information (Bulk Kit)


ARCIS DNA Sample Prep is a ready to use kit comprising two pre-filled tubes enabling pre-analytical processing of a variety of sample types including whole blood, saliva, bacteria, plant or animal cells or tissues.

In 3 minutes, with no prior sample preparation, the ARCIS DNA Sample Prep kit allows you to go from cells to downstream nucleic acid investigations without the need for isolation or purification.



2. Materials Provided

Material Provided	Quantity	Number of Preps
Tube 1: Lysis Buffer	1 Tube	48
Tube 2: Wash buffer	1 Tube	

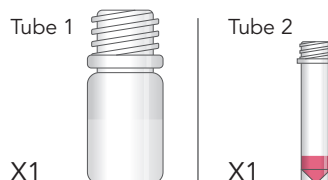

3. Storage Conditions

Tubes are shipped and stored at room temperature.

4. Applications

The product can be used on virtually any DNA containing sample. The nucleic acid released has been successfully applied to molecular biology techniques including PCR, sequencing and cloning.

The ARCIS DNA Sample Prep Kit accelerates pre-analytical processing in the following targets:

- Human Samples: including whole blood (fresh/frozen, with and without heparin or EDTA), plasma, urine and buccal swabs
- Microbiological samples: including bacteria (E.coli, S.aureus, P.aeruginosa, K. pneumoniae), viruses (HBV/HCV) and parasites (plasmodium)

5. Applications

The ARCIS DNA Sample Prep Kit is particularly suited to the following areas:

- Laboratory-free situations including point of care and field-based diagnostics, near-patient testing, and integrated sample-processing workflows.
- Forensic Analysis
- Veterinary research
- Genomics and biomarker analysis

Instructions for Use continued

6. Typical Protocol- Rapid processing of whole blood

Ensure samples have thawed completely before starting this procedure.

- 6.1. Add 30ul of whole blood to 150ul of Reagent 1 (or scale up for larger sample volume).
- 6.2. Incubate for one minute at room temperature.
- 6.3. Take 10ul of the above mixture and combine with 40ul of Reagent 2 (or scale up for larger sample volume).
- 6.4. Add appropriate volume into PCR master mix (e.g. 5µl per 25µl reaction).

For applications where enhanced sensitivity is required, the following protocol modifications can be used: Samples can be mixed with Reagent 1 at a 1:4 or 1:3 ratio to reduce sample dilution (See Table 1). Samples that have been processed in step 6.1 can be added to Reagent 2 at 1:3, 1:2 or 1:1 ratio to reduce sample dilution (See Table 2).

Table 1: Processing samples in Reagent 1

Sample Volume (µl)	Reagent 1 Volume (µl)	Ratio
30	120	1:4
60	240	1:4
90	360	1:4
30	90	1:3
60	180	1:3
90	270	1:3

Table 2: Washing samples in Reagent 2

Extract from Tube 1 (µl)	Reagent 2 Volume (µl)	Ratio
10	30	1:3
10	20	1:2
10	10	1:1

This product is for laboratory research use only.

CAUTION: Not for diagnostic use.

Dispose of any unused product within 48 hours after opening.

7. Manufacturer Contact Details

Arcis Biotechnology Ltd

The Innovation Centre
Sci-tech Daresbury
Keckwick Lane
Daresbury
Cheshire UK WA4 4FS
www.arcisbio.com



JH BIO Innovations Pvt. Ltd.

Ph: 080 - 23418944/45/46 Fax: 080- 23418947
E-Mail: info@jhindia.com Web: www.jhindia.com

